

LABORATORY-SCALE FLASH X-RAY DIFFRACTION

FROM SEEING MOTION
TO UNDERSTANDING
STRUCTURAL CHANGE,
NANOSECOND FLASH XRD
WITHOUT RELIANCE ON
LARGE-SCALE FACILITIES.

Flash XRD advantages


- Ultrafast pulse-resolved diffraction for dynamic processes
- Operando diffraction without pausing the experiment
- Direct access to phase, strain, and lattice evolution
- High-energy penetration in realistic sample geometries
- Precise timing and trigger-synchronized measurements
- Laboratory-scale alternative to synchrotron facilities

Example applications


Phase transition kinetics
Detection of combustion and chemical reaction
X-ray diffraction as a thermodynamic diagnostic
Dynamic compression studies
Rapid heating experiments
High strain-rate deformation analysis

Core components (fig 04)


Flash X-ray source
Electrodes
X-ray optics & collimator
Detector
Shielding and safety enclosure

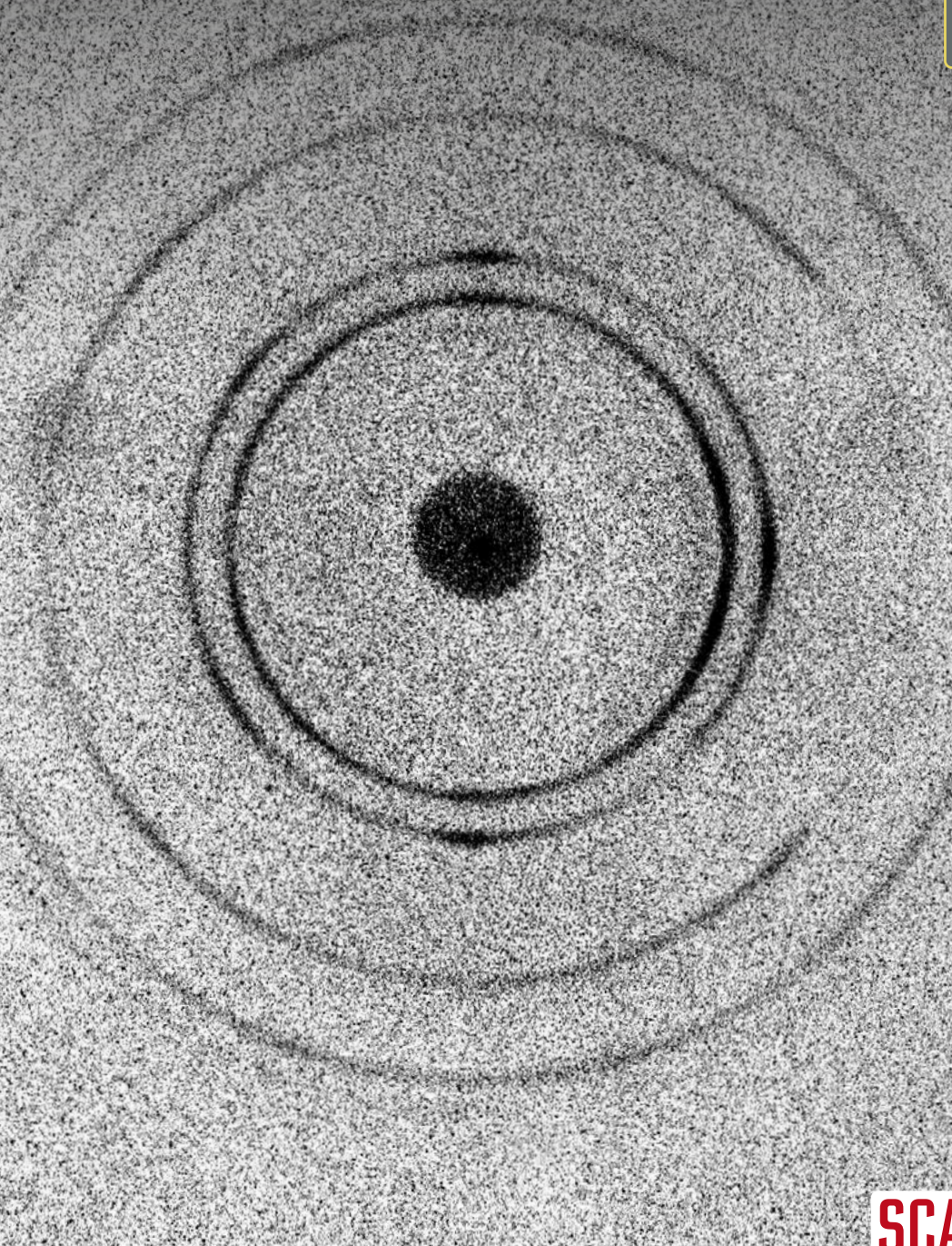


www.scandiflash.com

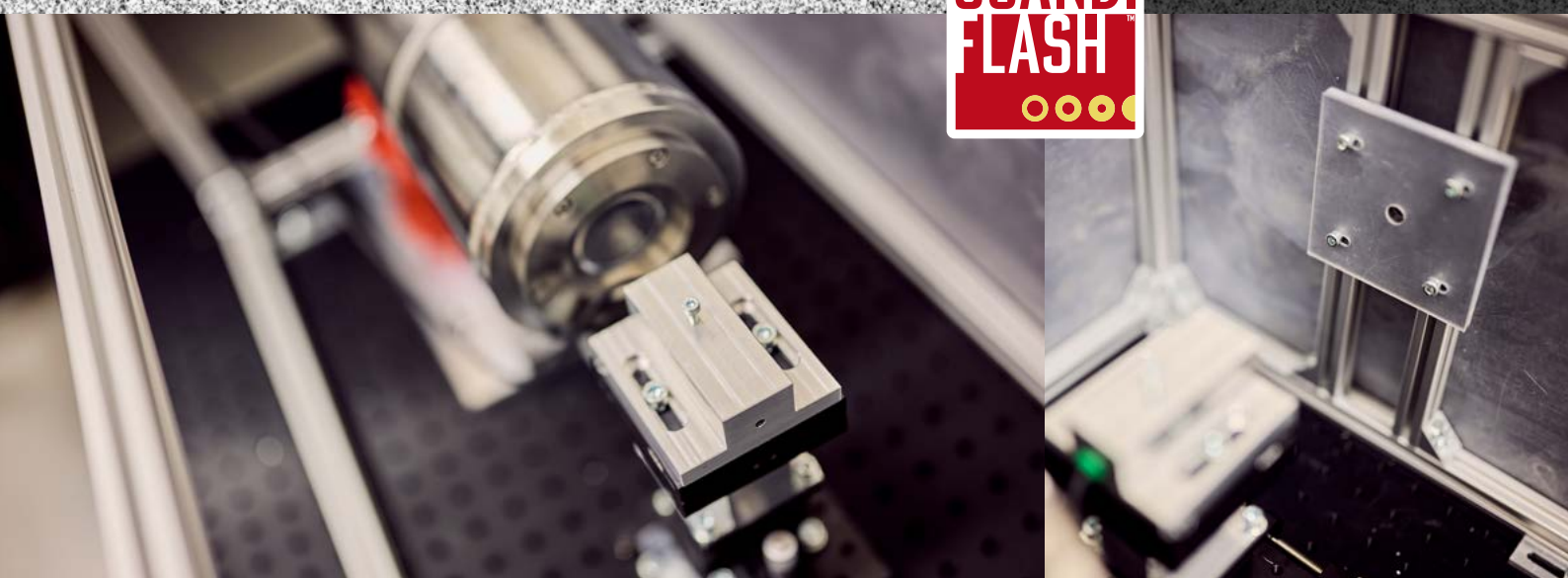


Raw XRD Curve (fig 01)

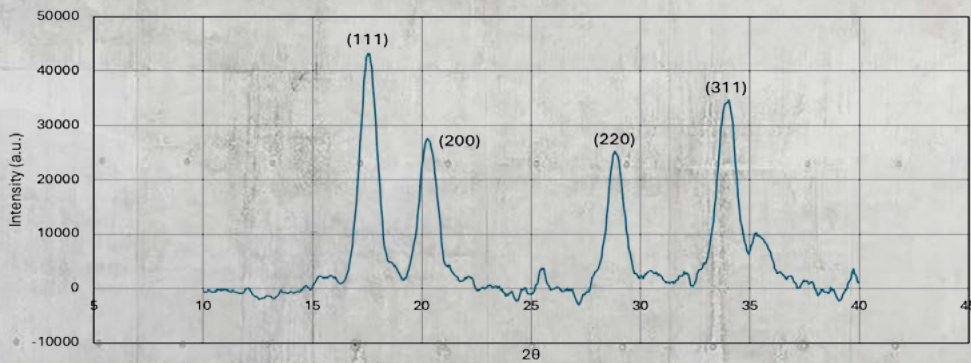


(fig 02)

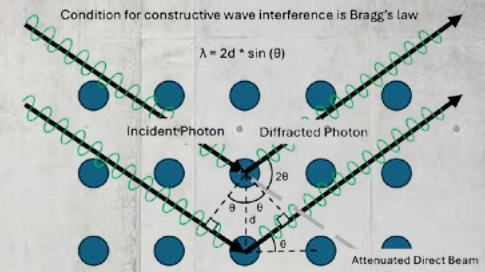

(fig 03)

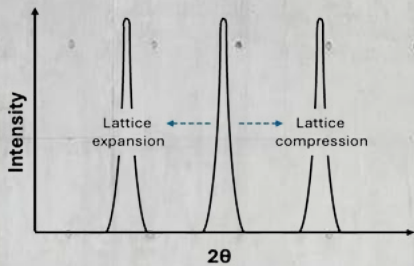

Component Setup (fig 04)

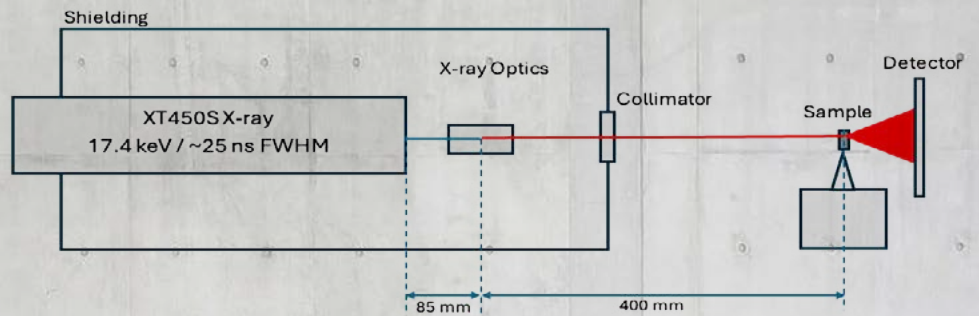

Understanding Structural Change in Real Time

X-ray diffraction (XRD) is one of the most widely used methods for studying crystalline materials. Because diffraction probes atomic lattice structure directly, it provides quantitative information about phase composition, lattice strain, and crystallographic orientation (fig 01).

While conventional laboratory XRD systems operate under steady-state conditions, flash X-ray diffraction enables measurements during transient events occurring on nanosecond to microsecond timescales. This makes Flash XRD a powerful diagnostic for dynamic materials research including shock compression, rapid heating, phase transitions, and high strain-rate deformation studies.

Structural Diagnostics During Dynamic Events

In a diffraction experiment, an incident X-ray beam interacts with lattice planes inside a crystalline material. When the Bragg condition is satisfied (fig 02), diffraction peaks or rings are produced that correspond to specific crystallographic planes.

Changes in diffraction peak position reveal changes in lattice spacing caused by compression, thermal expansion, elastic strain, and structural phase transitions. As a result, Flash XRD provides direct insight into the evolving thermodynamic and mechanical state of materials during dynamic experiments.

Time-Resolved Crystallographic Insight

Flash X-ray diffraction (Flash XRD) combines short, intense X-ray pulses with diffraction geometry to resolve structural evolution during fast, non-equilibrium processes. Unlike conventional laboratory XRD, which often requires long integration times and interrupted experiments, Flash XRD enables measurements while the process remains fully active — preserving the true material response.

This allows researchers to:

- Capture structural evolution in real time
- Reveal transient and intermediate states
- Understand cause-and-effect relationships
- Study ultrafast material processes
- Validate theoretical and computational models

From Shock Physics to Operando Materials Research

Flash XRD is well suited for shock physics, detonation studies, impact research, and other extreme-condition applications where structural changes occur on nanosecond timescales.

The technique also extends naturally to slower time-critical systems such as batteries and electrochemical devices, where the most important structural changes often occur during short transient events within otherwise long experimental cycles. By enabling operando diffraction during continuous operation, Flash XRD captures processes that conventional laboratory techniques may miss entirely.

Bridging the Gap Between Laboratory and Synchrotron XRD

Flash XRD bridges the gap between conventional laboratory diffraction and large-scale synchrotron experiments by combining crystallographic specificity with dramatically improved temporal resolution. The result is a compact laboratory-scale platform for dynamic diffraction studies that supports rapid experimental iteration, early-stage material screening, and synchronized multi-diagnostic research.

In many ways, Flash XRD is the structural analogue of high-speed imaging — not simply providing faster measurements, but enabling entirely new classes of observable material behavior.

Diffraction as a Thermodynamic Diagnostic

X-ray diffraction provides a non-contact diagnostic capable of inferring thermodynamic variables through measurements of lattice spacing.

Compression shifts diffraction peaks toward higher 2θ angles, while heating and expansion shift peaks toward lower angles (fig 03).

These peak shifts are governed by the material equation of state (EOS), enabling diffraction measurements to diagnose pressure, temperature, strain, and density evolution during dynamic experiments.



scan this for more info